



Profilometer User Operation manual

1. Log in on the profilometer computer by using :
User name: .\Microfluidics User
Password: Goldenson117
2. Turn on the Profilometer: On/off switch located on the back of the equipment:

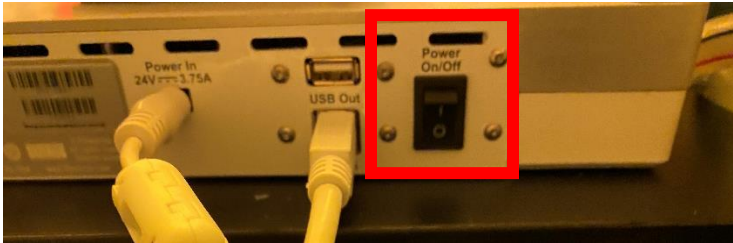


3. Open profilometer application: “**Profilm**”  located on the desktop or by typing on the search bar.


4. Once the program is open, wait for the auto calibration to see the Live window:

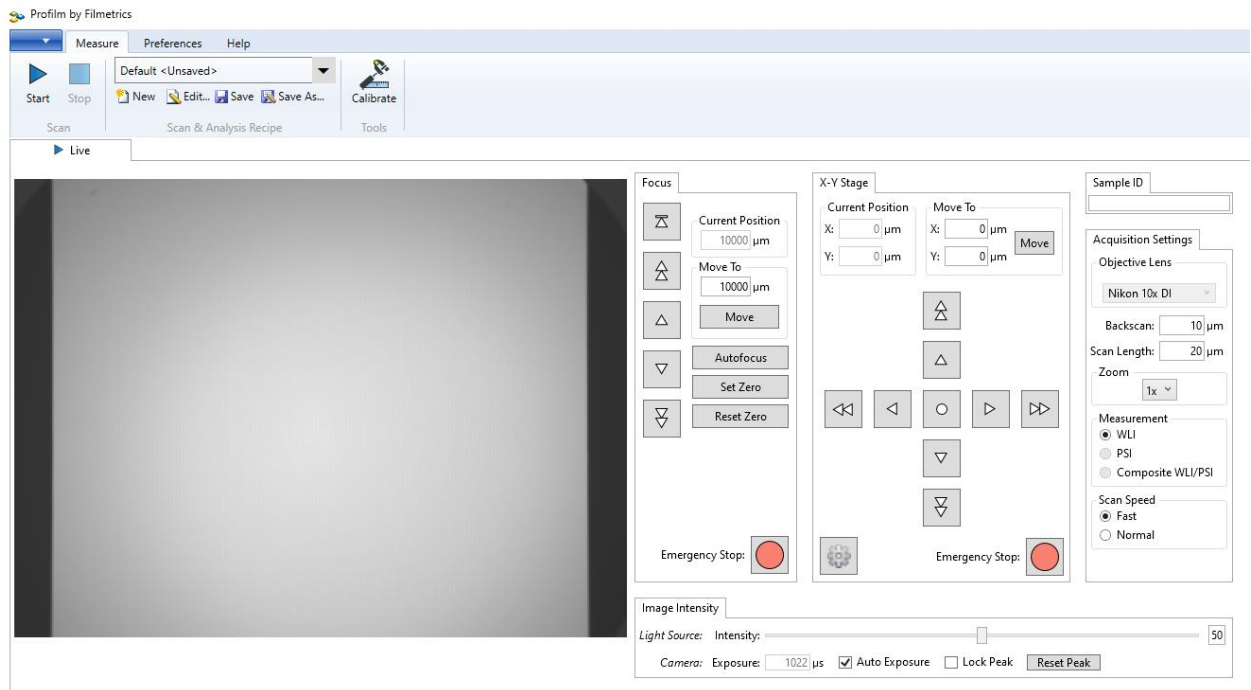


- Place the Master/PDMS that will be measured on the main stage. Adjust the controls (focus up and down and platform current motion X,Y) on the live window to have the desire image.

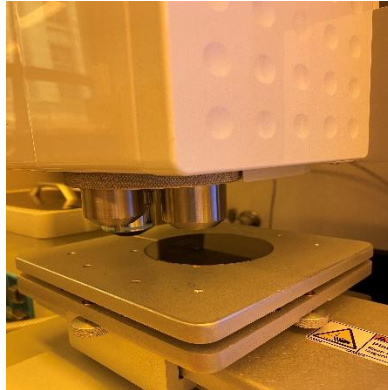


- Once desire image is achieved, click on the Start button in the live window to start the scan:

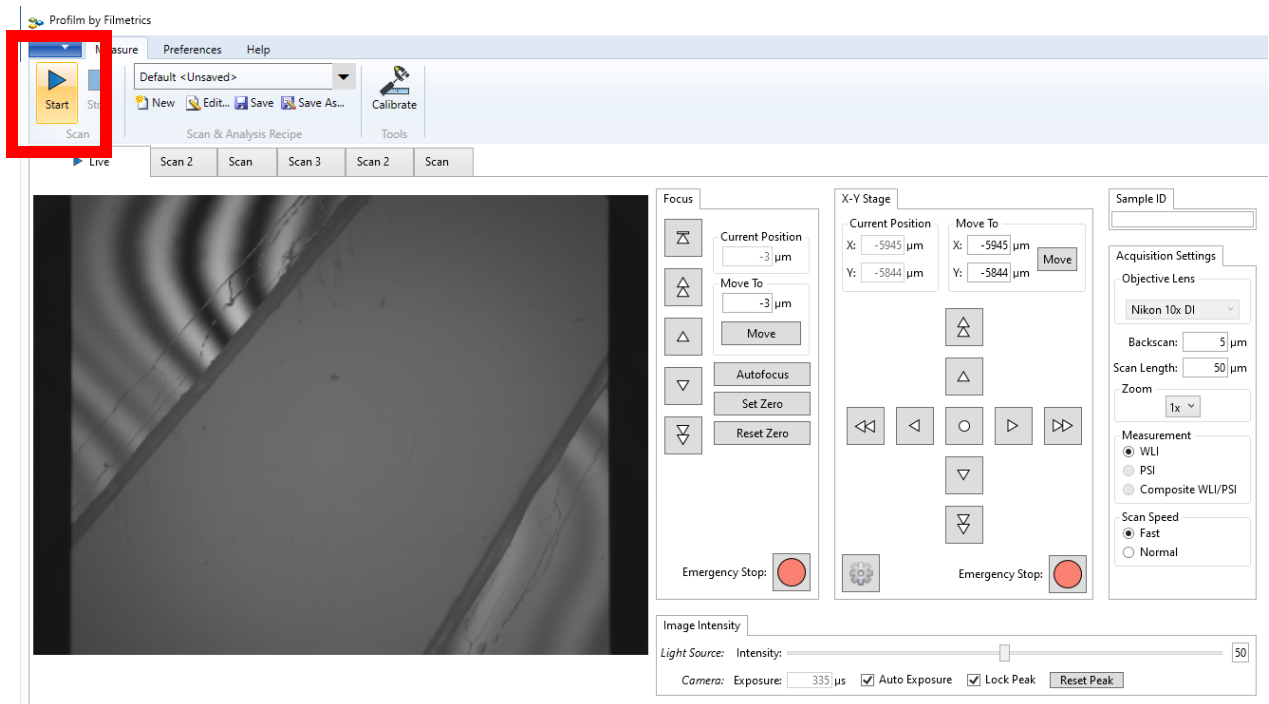


7. Once a scan is completed, it will open as an individual tab:

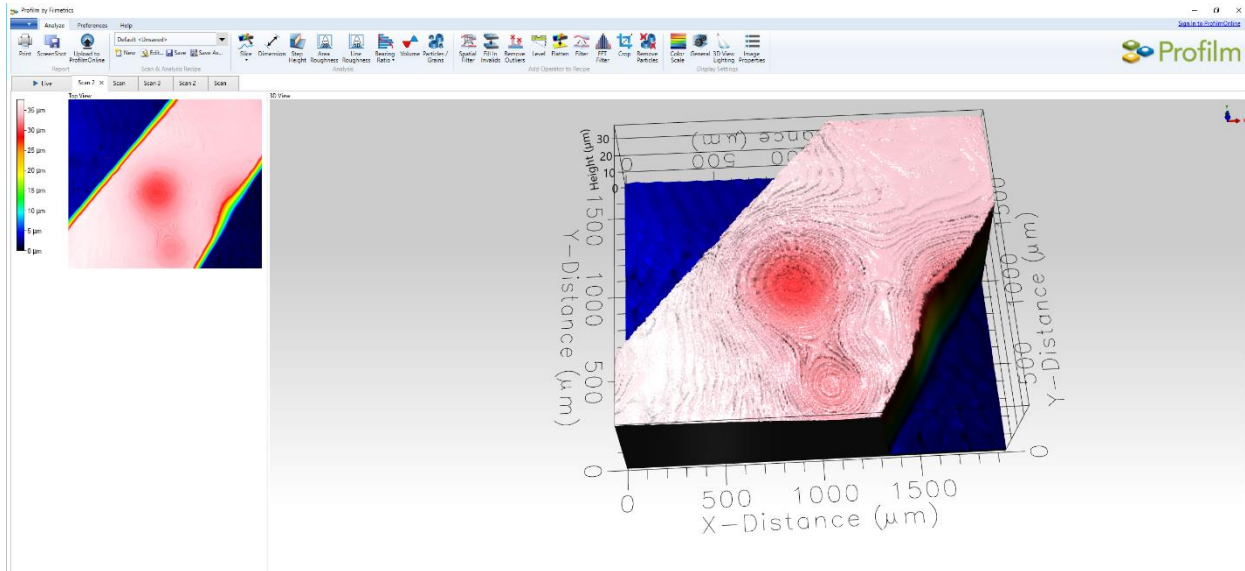

8. After the desired scan is obtained, you can use any of the controls/features located on the top of the image window for analysis:

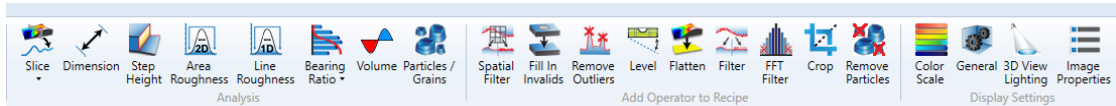

9. To start a new scan, go back to the live window and repeat the process.

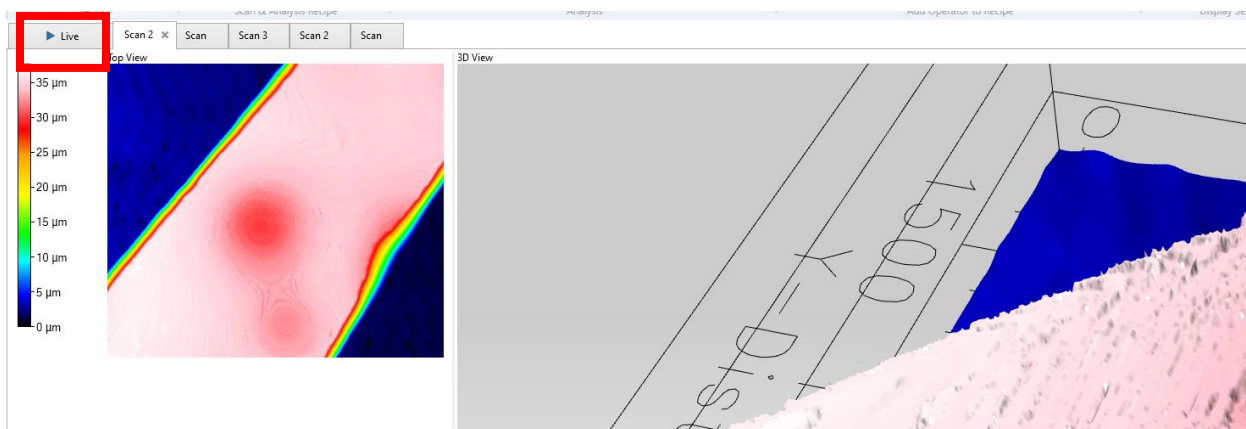

10. Once done using the application, close the program.

11. Turn off the Profilometer: On/off switch located on the back of the equipment.

12. Log out from the computer.